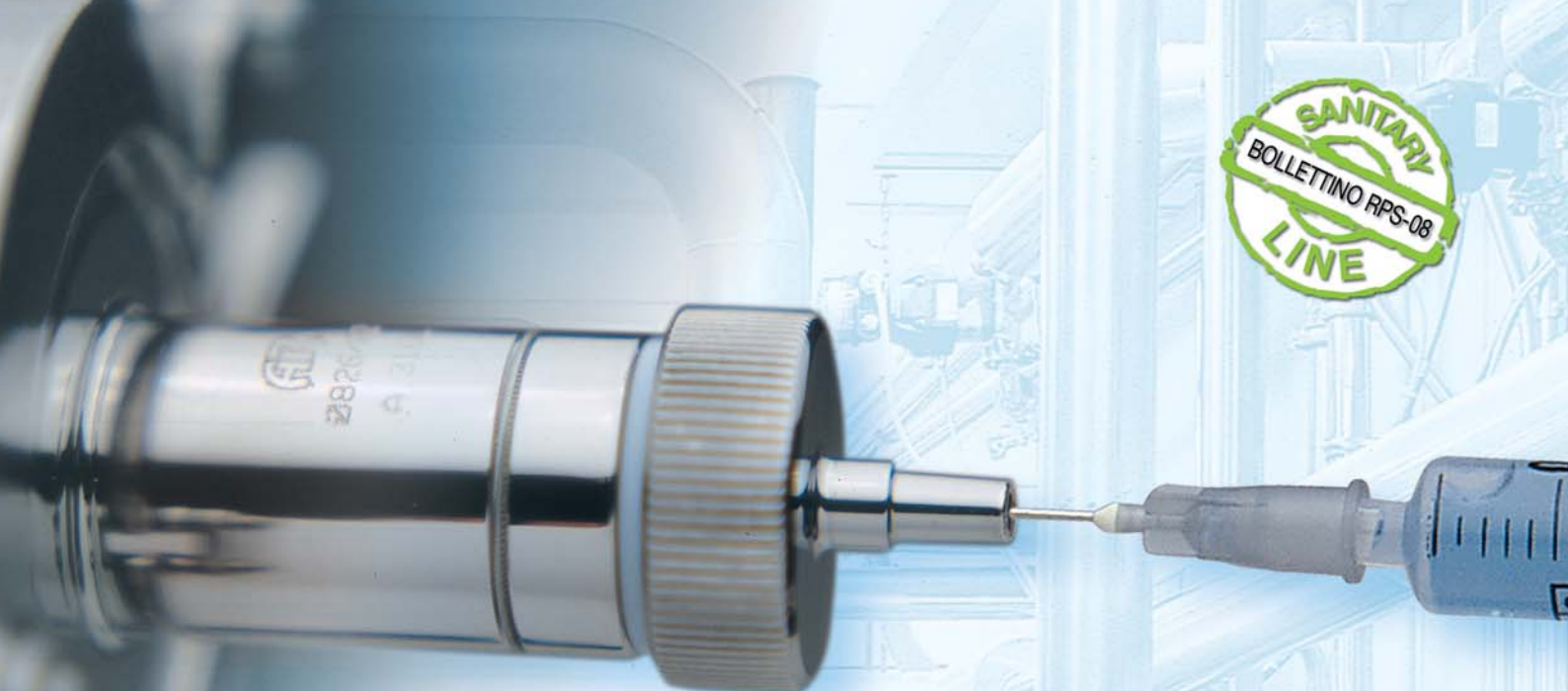
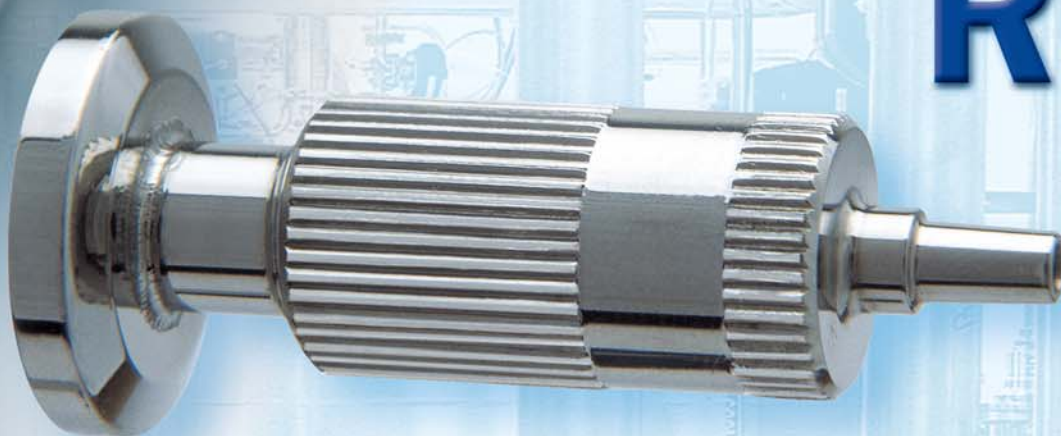
RPS

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**SANITARY
SAMPLING
VALVES**

Rps Available Body

**Standard
CLAMP 3/4" BS**

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RPS SAMPLING

CHARACTERISTICS

The **RPS** fully autoclavable sampling valves are designed for aseptic sampling of liquids and pharmaceutical products.

Made entirely in AISI 316L, the standard version is in the form of a rod with a 3/4" clamp and no welding.

On request, the connection to the process may be modified as indicated.

Easily installed onto distribution lines or onto tanks, this valve can be kept in aseptic conditions between one sample and another by using a silicon cap which closes the "luer slip" outlet. It's special design prevents deposits of bacteria and particles, possibly distorting the results of bacteriological tests carried out while allowing samples to be taken from the middle of the line enabling a representative sample to be taken. Extremely economical and very easy to use, the valve opens by turning the head knurled ratchet in an anticlockwise direction and then clockwise to return it to the closed position.

TECHNICAL DATA

MATERIALS

Valve body: AISI 316L

Shaft: AISI 316L

O-ring: Standard silicone
Optional viton

Plug: Standard silicone
Optional viton or PTFE

MAX WORKING CONDITIONS

Max Pressure: to 10 bar and 150°C

PROCESS CONNECTIONS

Standard: clamp 3/4" BS

Optional: clamp 1" BS
3/8"-1/2" gas male thread
butt-weld

OUTLET CONNECTIONS

Standard: LUER SLIP joint

Optional: Rubber hose Ø 6mm
Rapid connection

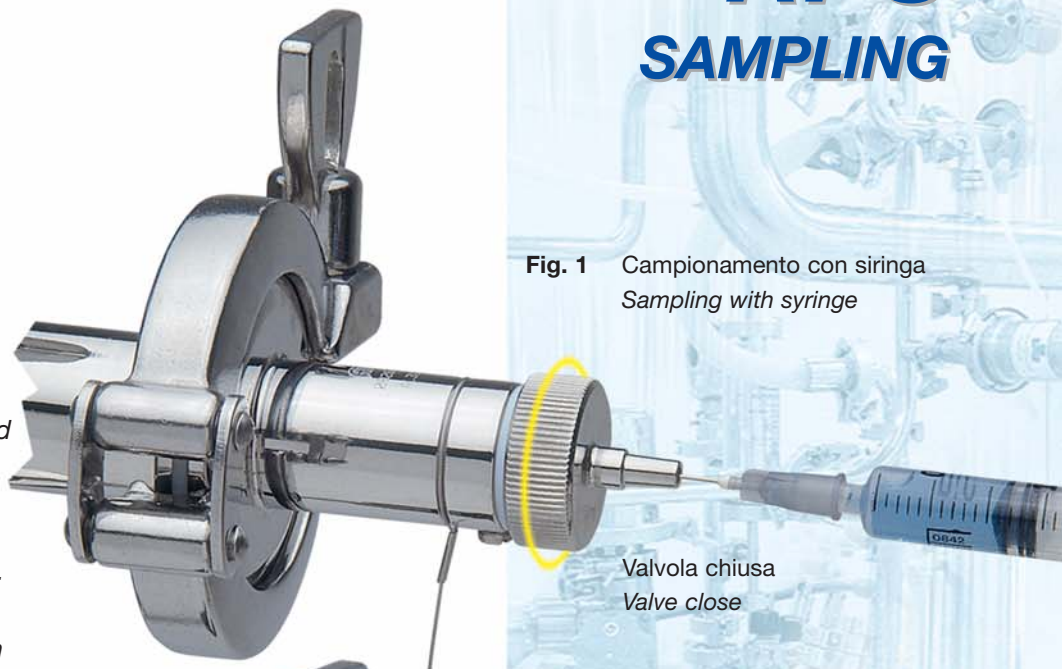
Fig. 1 Campionamento con siringa
Sampling with syringe

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
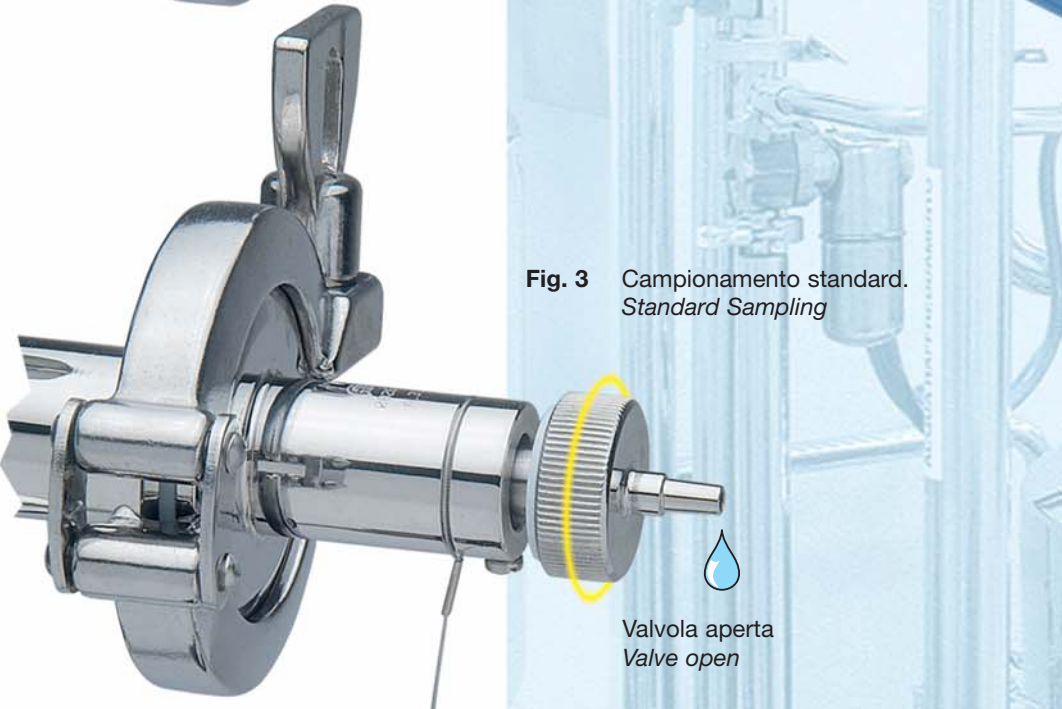
Fig. 3 Campionamento standard.
Standard Sampling

Valvola aperta
Valve open

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
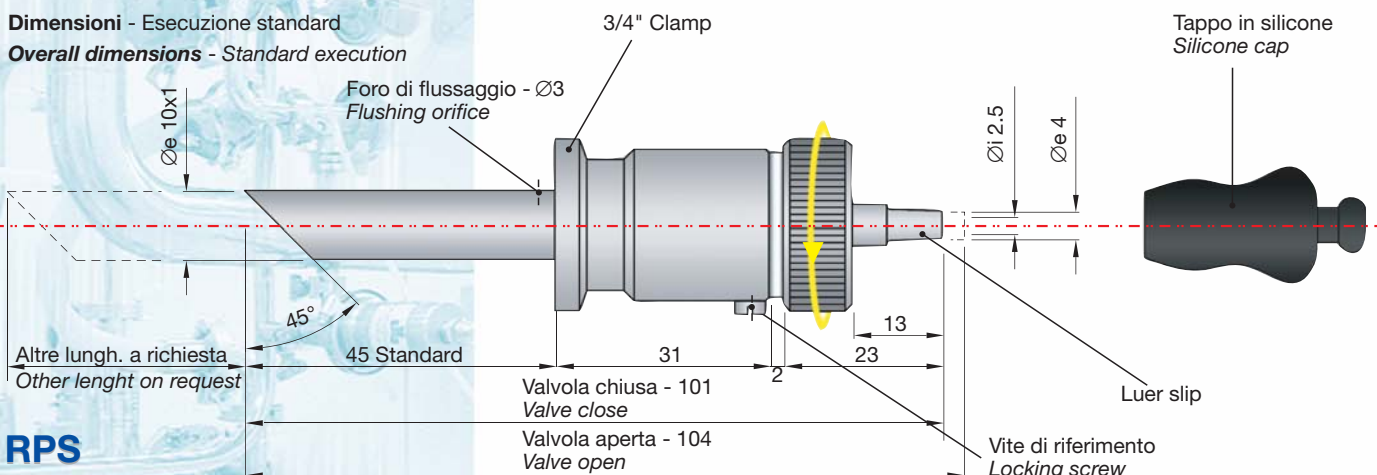
Overall dimensions - Standard execution

RPS

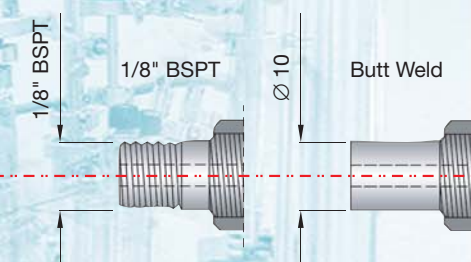RPS-VENT

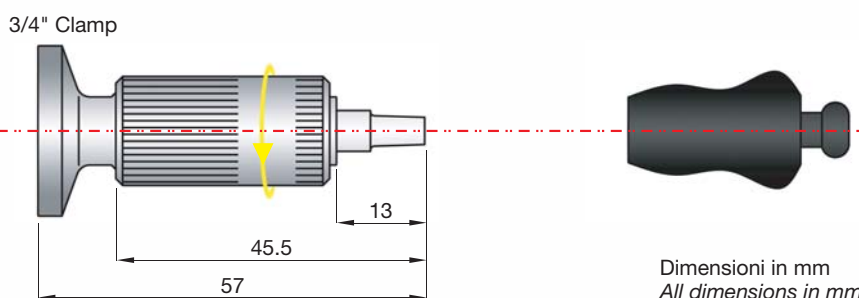Dimensioni in mm
All dimensions in mm

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INSTALLATION

Depending on the type of connection, the valve may be installed by welding, screwing or clamping the body. As the direction of flow is one-way, check the direction of the reference screws to ensure the body valve flow is correct. For all types of connection, ensure the luer outlet is horizontal to the valve and parallel with the floor. If flow is vertically upwards, the reference screws must be placed at the bottom, while if the flow is vertically downwards, the reference screws must be placed at the top.

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ATTACCO RAPIDO - optional
RAPID CONNECTION - optional

PORTAGOMMA - optional
RUBBER HOSE - optional

Sanitary standards symbol
administrative council
certificate of authorization # 1037
for standard no. 59-00

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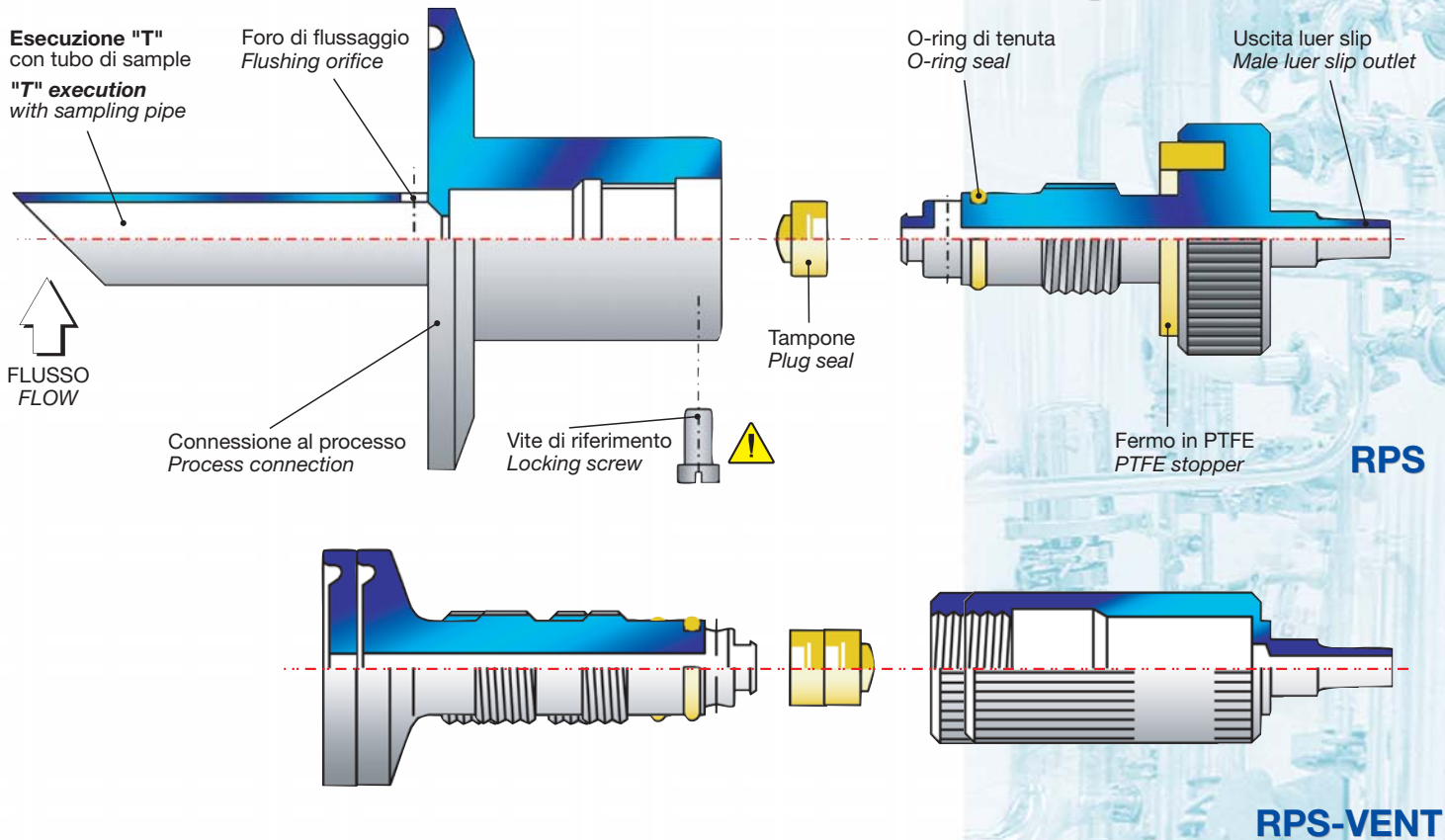
DOCUMENTATION

All valves are fully traceable for validation processes. Each valves is accompanied by material test certificates EN 10 204 3.1B 3A standards compliance certificates, FDA regulations compliance certificates, and tapes detailing the valve interior surfaces Ra.

Sanitary sampling valve

Ricambi Spare parts

cod. TRPS-V

Tampone e o-ring in VITON - siringabile
Plug seal and o-ring in VITON - syringable

cod. TRPS-X

Tampone e o-ring in SILICONE - siringabile
Plug seal and o-ring in SILICONE - syringable

cod. TRPS-G

Tampone in PTFE e o-ring in VITON - NON siringabile
Plug seal in PTFE and o-ring in VITON - NOT syringable

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H faccia piana
flat face

Attacco sample
Sample connection
LUER luer slip
P portagomma
rubber hose
S attacco rapido
rapid connection

LUER luer slip
P/S

VPA - VPAK
ASEPTIC SAMPLING VALVES

RPS
SANITARY SAMPLING VALVES

SSB
SANITARY SAMPLING BOTTLE

VA - VFA
TANK BOTTOM VALVES

VRN
SPRING CHECK VALVES

VSS
HIGH PURITY BALL VALVES

VF
BUTTERFLY VALVES

HE
DTS HEAT EXCHANGERS

SP - SL
SIGHT GLASS-FLOW INDICATOR

TC
CLAMP FITTINGS

TK - CONN
TANK CONNECTIONS

RE - FLEX
SILICONE HOSE & FITTINGS

MM
MAGNETIC MIXER

DD
REPLACEMENT DIAPHRAGMS

RSH
ROTATIVE SPRAY HEAD

**Sanitary
flow
equipment**