



Ezi-Flow CSV Chargebag filling Operating Instructions

Follow these instructions for manually filling
an Ezi-Flow CSV Chargebag



THE QUEEN'S AWARDS
FOR ENTERPRISE:
INNOVATION
2014



THE QUEEN'S AWARDS
FOR ENTERPRISE:
INNOVATION
2015



THE QUEEN'S AWARDS
FOR ENTERPRISE:
INNOVATION
2019

VERSION 1.0
DATE PRODUCED 04/03/22


1

- During this procedure
- you will need:

- 1 Ezi-Flow CSV Chargebag with a pre-fitted Passive
- 2 Ezi-Grip Single-Use Bag Clamp
- 3 Red Tamper Evident Security Tag
- 4 Filling Funnel – either Stainless Steel or Polypropylene
- 5 CSV Passive Opening Tool

In addition, a number of mobile or benchtop filling stands are also available to make the Filling Process easier.



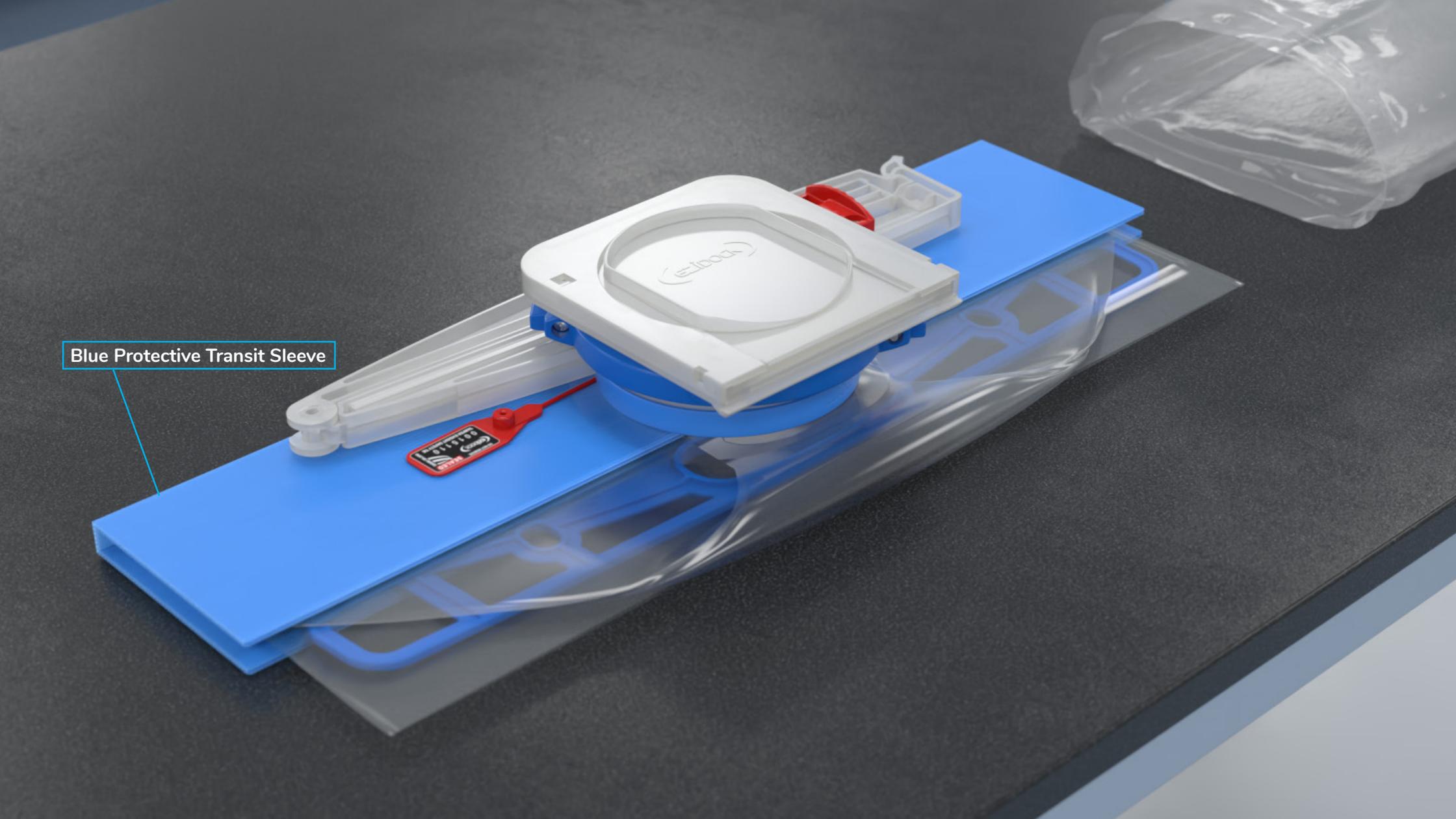


Blue Protective Transit Sleeve

2:

First, remove the Chargebag, Red Tamper Evident Security Tag and Ezi-Grip Single-Use Bag Clamp from the pouch by tearing the Perforated Seal. The Blue Protective Transit Sleeve can be discarded.

EZI-DOCK SYSTEMS LTD
Ezi-Flow CSV
Chargebag Filling

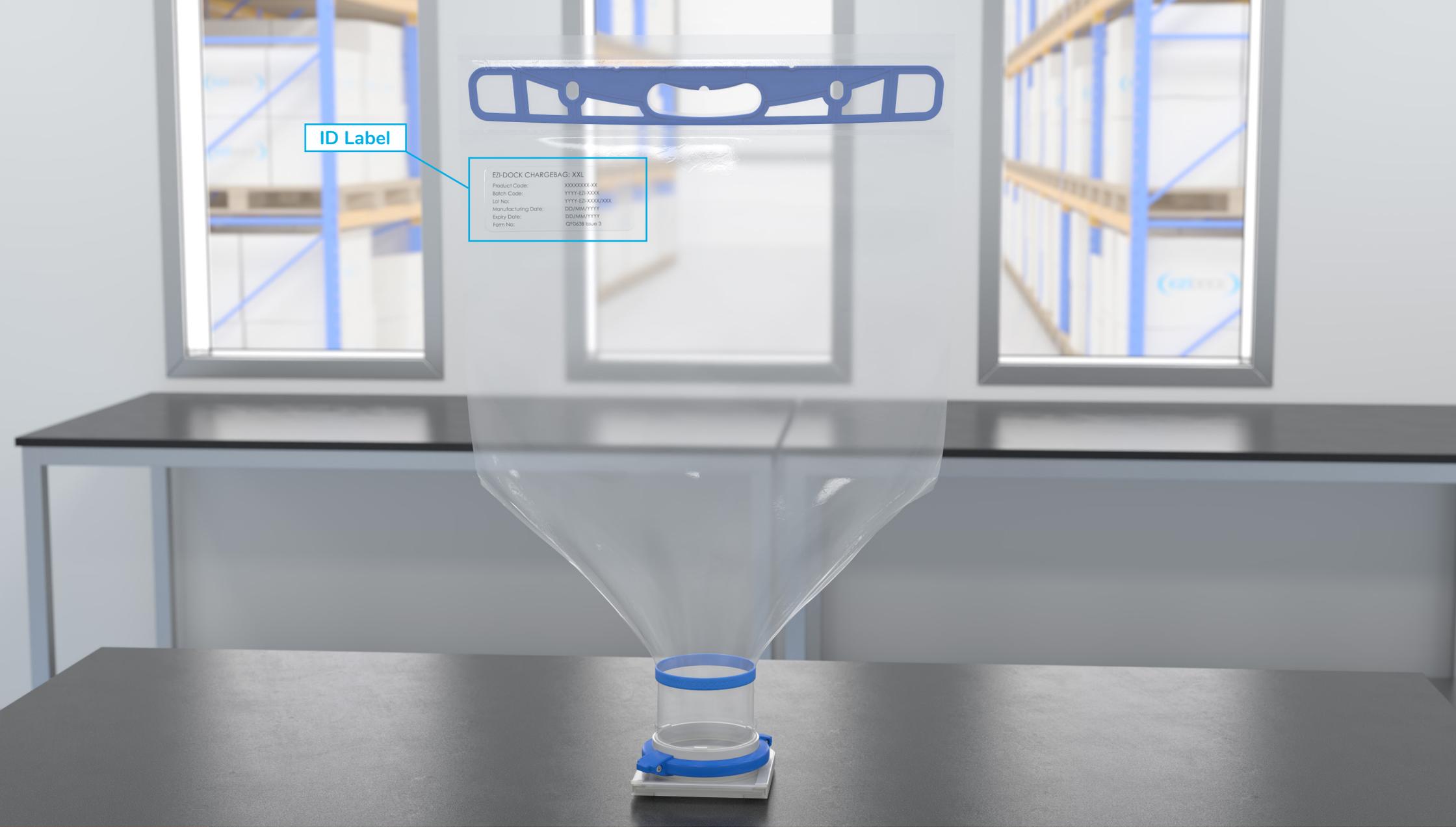


ID Label

EZI-DOCK CHARGEBAG: XXL
Product Code: XXXXXXXXXX
Batch Code: YYYY-EZ-XXXX
Lot No: YYYY-EZ-XXXX/XXX
Manufacturing Date: DD/MM/YYYY
Expiry Date: DD/MM/YYYY
Form No: GP0638 Issue 3

3:

Each Chargebag has a Clear ID Label showing all Manufacturing Details, including Batch & Lot Number, Part Number and Chargebag Volume, as well as Production and Expiry dates.

EZI-DOCK SYSTEMS LTD
Ezi-Flow CSV
Chargebag Filling

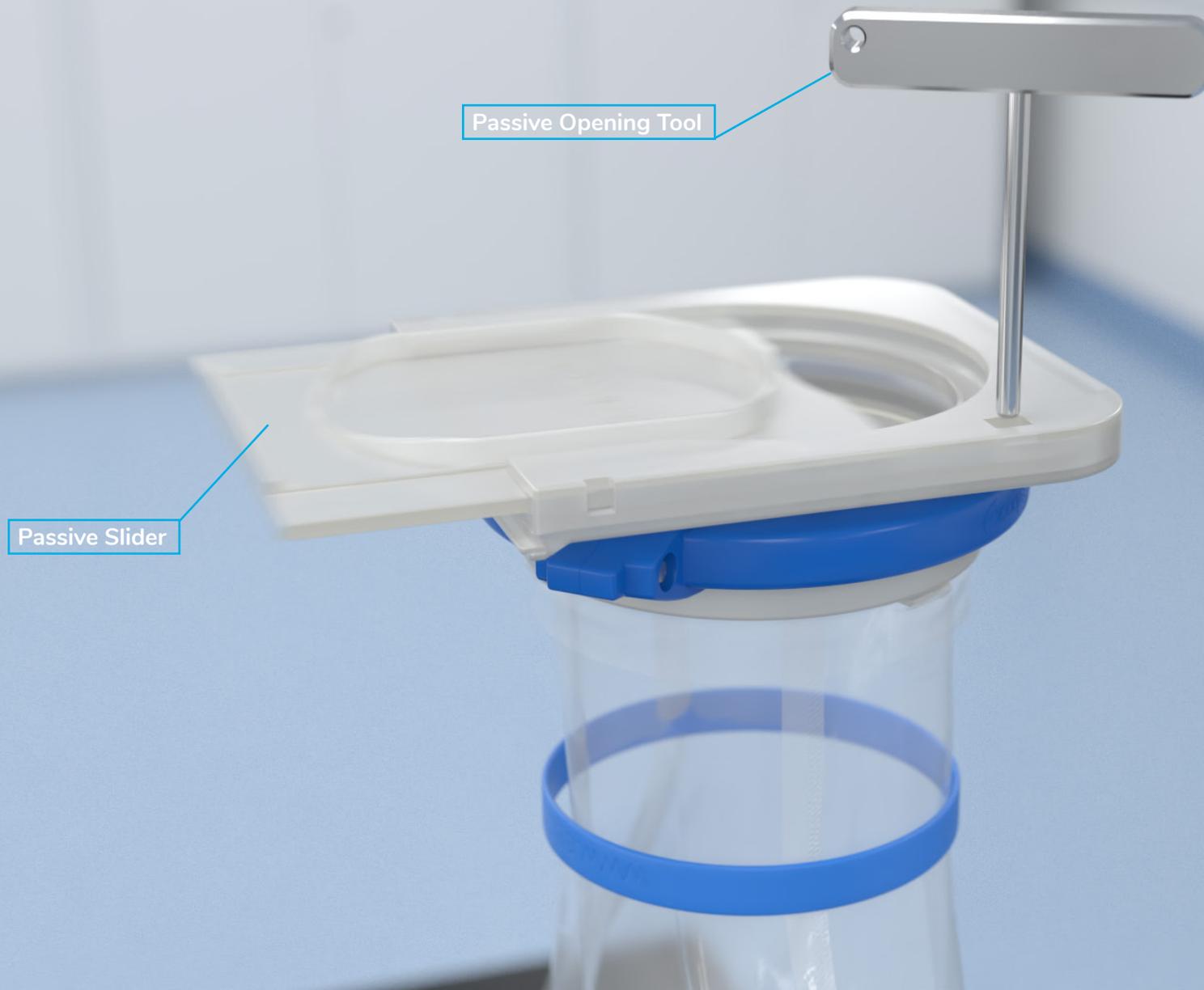

4:


The Ezi-Flow CSV Passive has a Manual Interlock which prevents the Passive Slider from being removed. To remove the Passive Slider, simply depress the Passive Interlock Mechanism using the Passive Opening Tool. Now you can fully withdraw the Slider from the passive Body.



Filling Stand

5:

Next, place the Chargebag into the Filling Stand with the ID Label facing forwards.

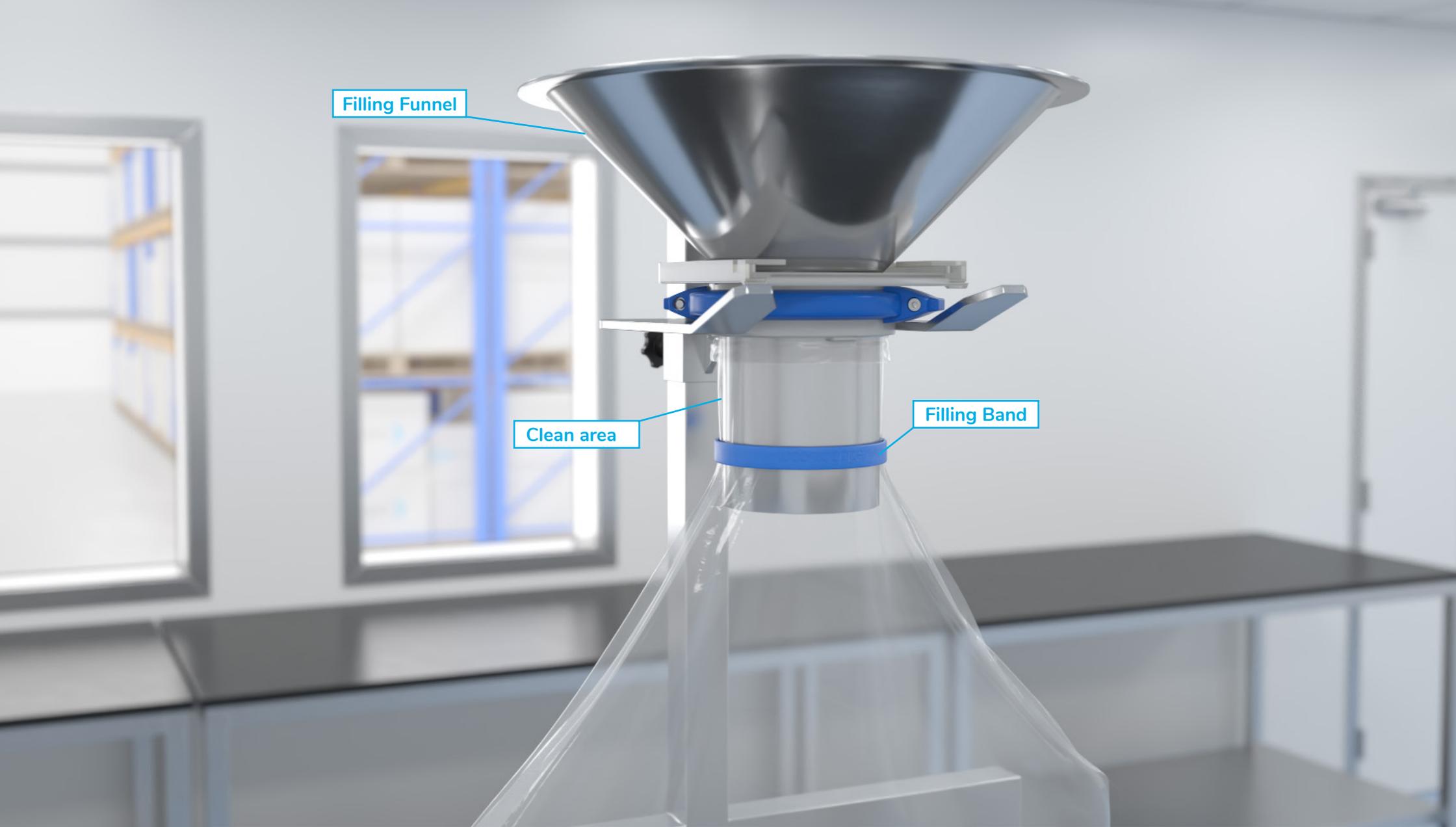

Filling Funnel

Clean area

Filling Band

6:

Fully insert the Filling Funnel through the Passive Body. The Filling Band must tightly wrap around the Filling Funnel and should be positioned towards the end of the Funnel Spout. This will create a “Clean Area” and prevent product from contaminating the Chargebag Neck & Passive during the Filling Process.

EZI-DOCK SYSTEMS LTD
Ezi-Flow CSV
Chargebag Filling



7:

Product can now be carefully dispensed through the Filling Funnel into the Chargebag.


EZI-DOCK SYSTEMS LTD
Ezi-Flow CSV
Chargebag Filling



8:

When Filling is complete, carefully remove the Funnel, ensuring that no residual product is retained on the surface of the Funnel.

EZI-DOCK SYSTEMS LTD
Ezi-Flow CSV
Chargebag Filling

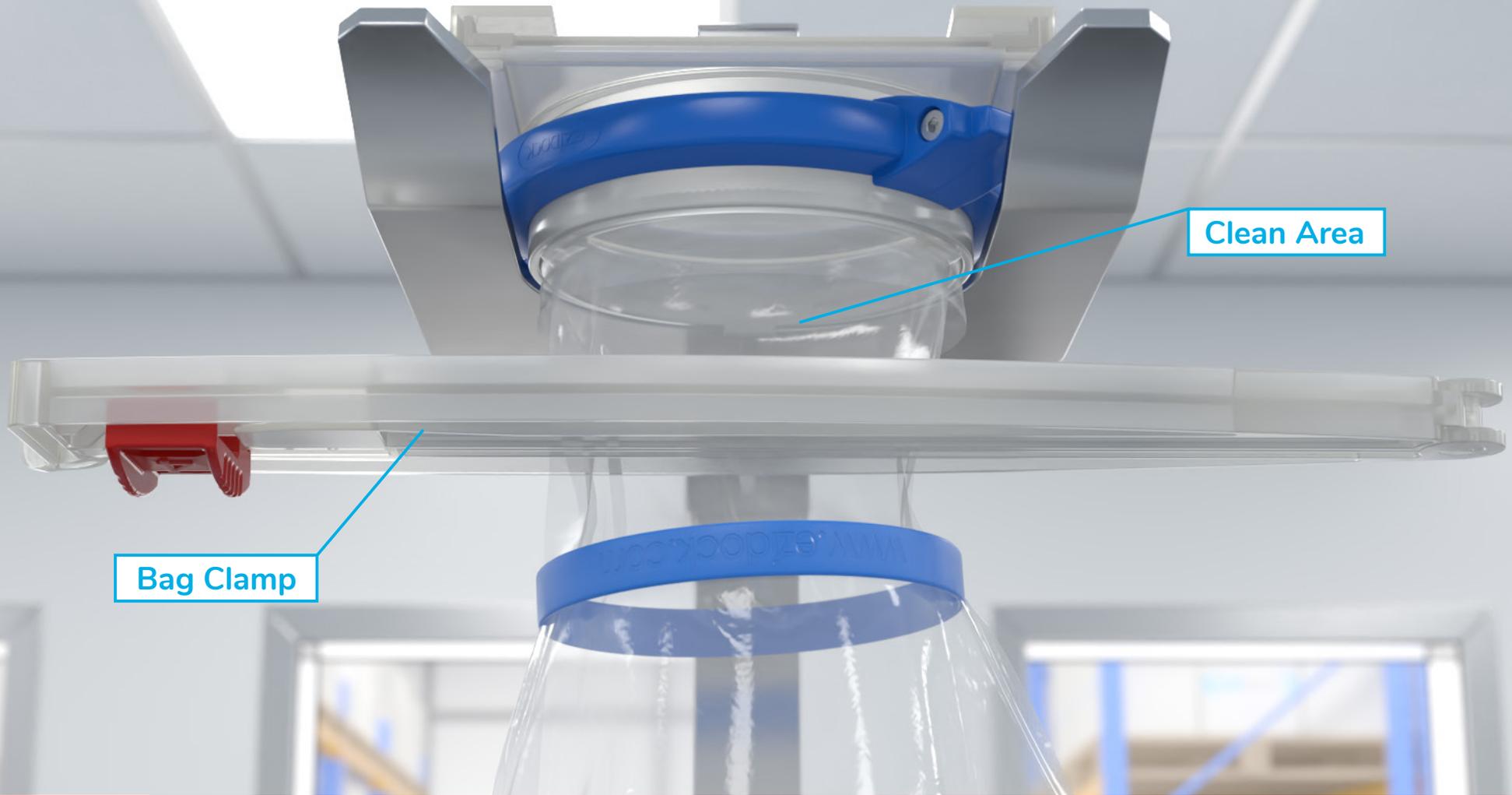


9:

Next, the Ezi-Grip Single-Use Bag Clamp needs to be used to seal off the neck of the Chargebag and prevent product from contaminating the Passive. Be sure to position the Clamp in the “Clean Area” between the Passive and the Filling Band.



Ensure the clamp is orientated with the red toggle facing downwards and to the left. This is an important step and best working practice when using the system.

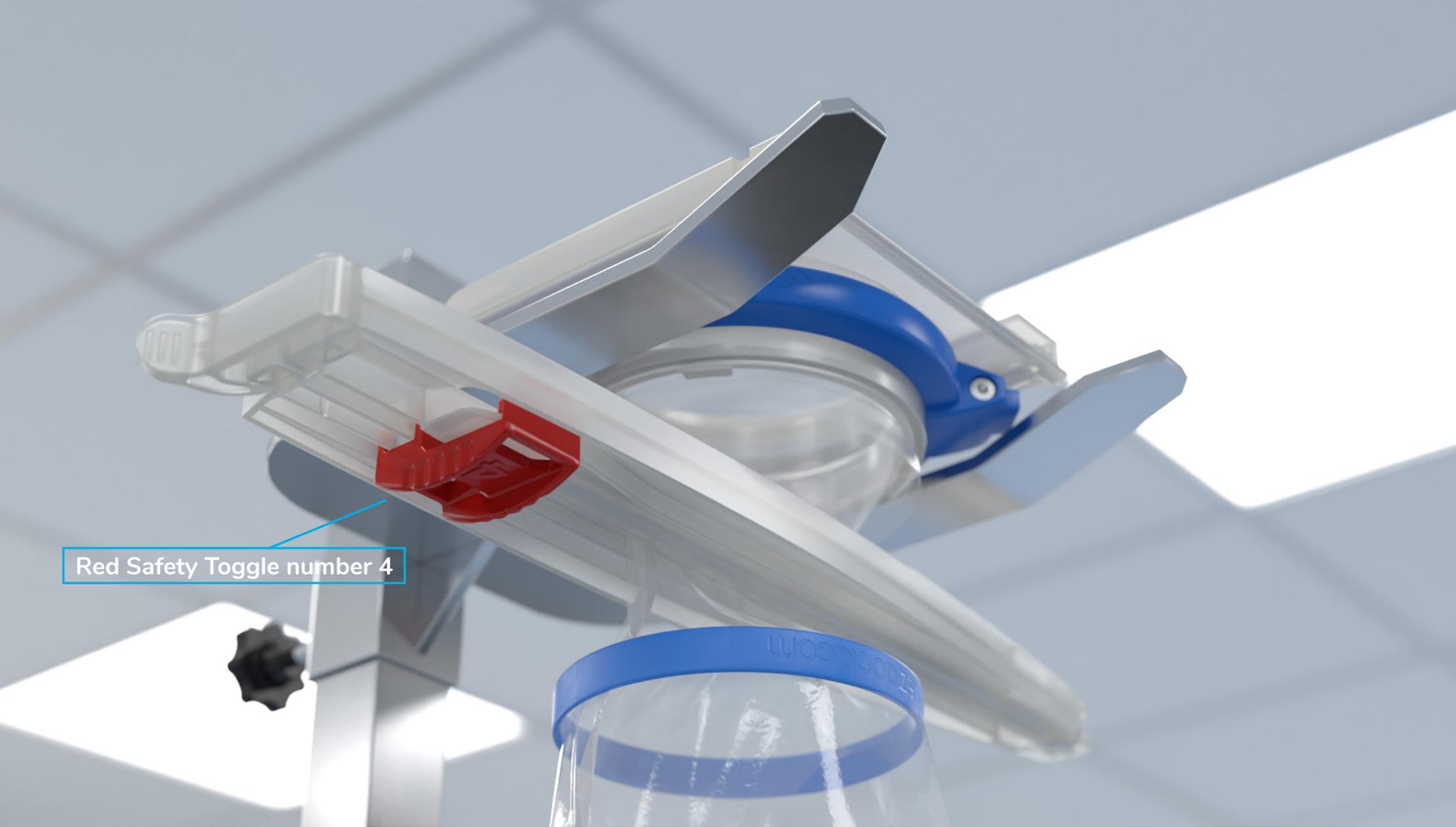


Red Safety Toggle number 4

10:

Carefully, push the two halves of the Bag Clamp together.

When they are fully engaged, slide the Red Safety Toggle, marked number 4, towards the centre of the clamp.

EZI-DOCK SYSTEMS LTD
Ezi-Flow CSV
Chargebag Filling

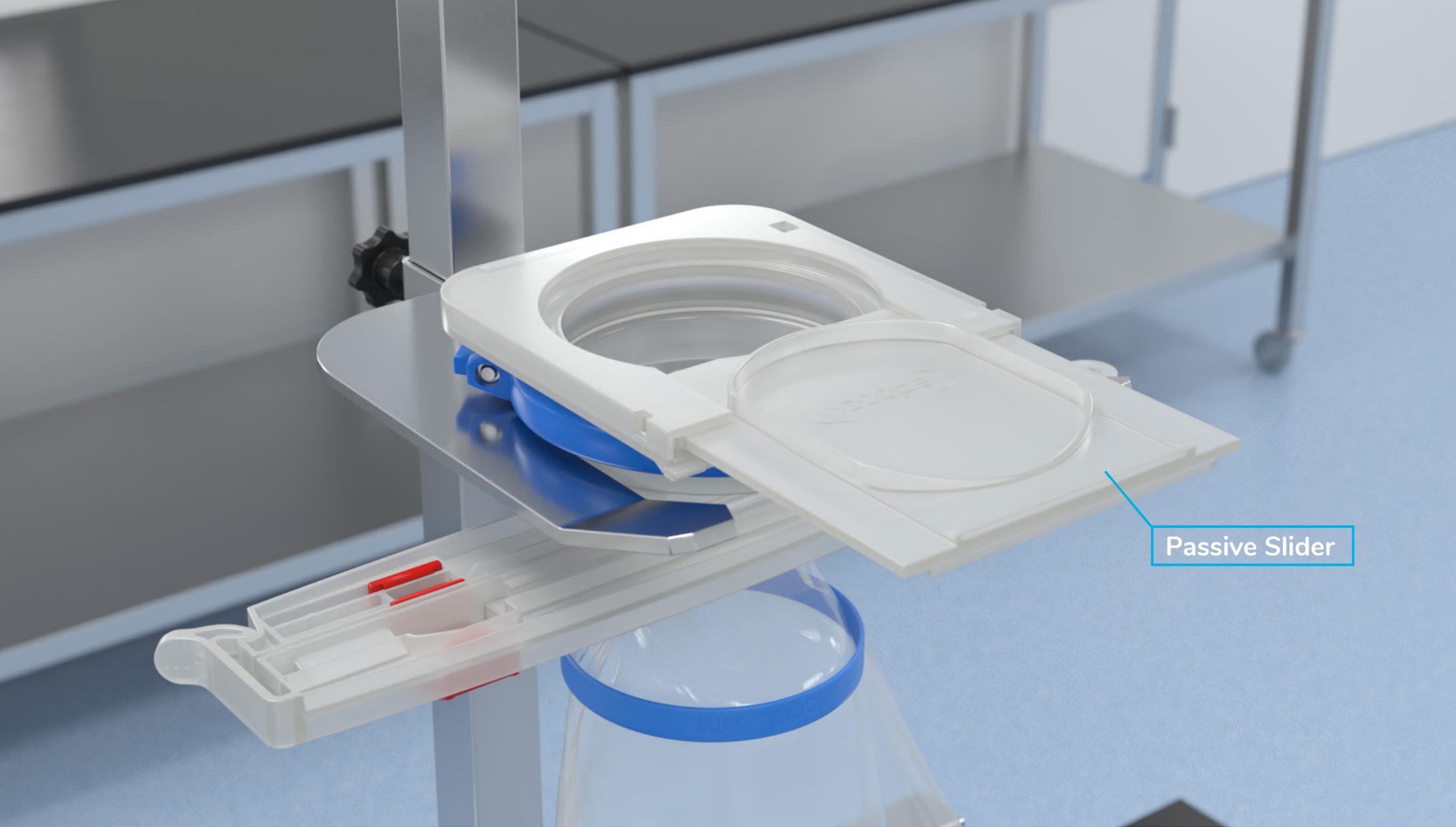


Passive Slider

11:

The Passive Slider can now be re-inserted into the Passive Body. This will re-engage the Interlock.

EZI-DOCK SYSTEMS LTD
Ezi-Flow CSV
Chargebag Filling

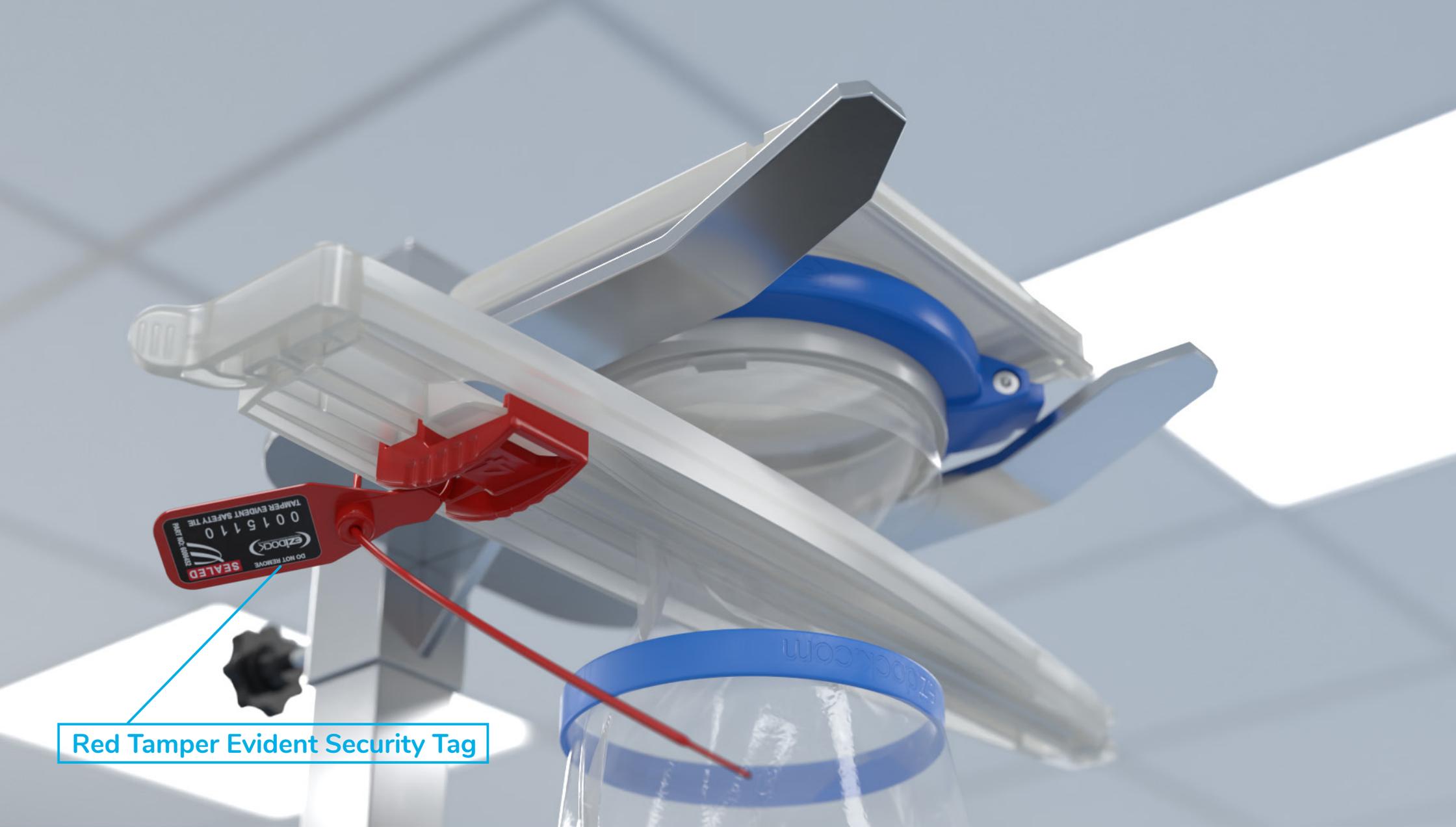

Red Tamper Evident Security Tag

12:

Finally, insert the Tamper Evident Security Tag into the holes in the Clamp Handle and pull tightly. This will prevent accidental release of the Bag Clamp.

EZI-DOCK SYSTEMS LTD
Ezi-Flow CSV
Chargebag Filling



13:

The filling procedure is now complete, and the Chargebag is ready for discharge.

EZI-DOCK SYSTEMS LTD
Ezi-Flow CSV
Chargebag Filling



THE QUEEN'S AWARDS
FOR ENTERPRISE:
INNOVATION
2014



THE QUEEN'S AWARDS
FOR ENTERPRISE:
INNOVATION
2015



THE QUEEN'S AWARDS
FOR ENTERPRISE:
INNOVATION
2019

Ezi-Dock Systems Limited

Park Court, Park Lane Business Park, Kirkby in Ashfield,
Notts NG17 9GU United Kingdom

 Tel: **+ 44 (0) 1623 888 000**

 Email: **sales@ezidock.com**

 Web: **www.ezidock.com**